

SuperMOUNT™ Mounting Medium

CATALOG #: 1211-20

AMOUNT: 20 ml

LOT #: _____

STORAGE CONDITIONS: Store at -20°C.

SHELF LIFE: Stable for up to 6 months at -20°C

DESCRIPTION:

The SuperMount™ Mounting Medium was prepared using a unique, stable formula to preserving fluorescence of fluorescein, rhodamine, Texas Red as well as other fluorescent reagents. Sections mounted in the Mounting Medium retain fluorescence during prolonged storage.

RECOMMENDED USAGE:

To mount tissues or cells on a slide, dispense one drop of the Mounting Medium onto the section. Then coverslip and allow the mounting medium to disperse over the entire section. The SuperMount Mounting Medium does not solidify, but remains a liquid on the slide. After mounting, coverslipped slides will not readily dry out.

For prolonged storage coverslips can be permanently sealed around the perimeter with nail polish or a plastic sealant. Mounted slides should be stored at 4°C protected from light.

FOR RESEARCH USE ONLY! Not to be used in humans.

RELATED PRODUCTS:

Apoptosis Detection Kits & Reagents

- Annexin V Kits & Bulk Reagents
- Caspase Assay Kits & Reagents
- Mitochondrial Apoptosis Kits & Reagents
- Nuclear Apoptosis Kits & Reagents
- Apoptosis Inducers and Set
- Apoptosis siRNA Vectors

Cell Fractionation System

- Mitochondria/Cytosol Fractionation Kit
- Nuclear/Cytosol Fractionation Kit
- Membrane Protein Extraction Kit
- Cytosol/Particulate Rapid Separation Kit
- Mammalian Cell Extraction Kit
- FractionPREP Fractionation System

Cell Proliferation & Senescence

- Quick Cell Proliferation Assay Kit
- Senescence Detection Kit
- High Throughput Apoptosis/Cell Viability Assay Kits
- LDH-Cytotoxicity Assay Kit
- Bioluminescence Cytotoxicity Assay Kit
- Live/Dead Cell Staining Kit

Cell Damage & Repair

- HDAC Fluorometric & Colorimetric Assays & Drug Discovery Kits
- HAT Colorimetric Assay Kit & Reagents
- DNA Damage Quantification Kit
- Glutathione & Nitric Oxide Fluorometric & Colorimetric Assay Kits

Signal Transduction

- cAMP & cGMP Assay Kits
- Akt & JNK Activity Assay Kits
- Beta-Secretase Activity Assay Kit

Adipocyte & Lipid Transfer

- Recombinant Adiponectin, Survivin, & Leptin
- CETP Activity Assay & Drug Discovery Kits
- PLTP Activity Assay & Drug Discovery Kits
- Total Cholesterol Quantification Kit

Molecular Biology & Reporter Assays

- siRNA Vectors
- Cloning Insert Quick Screening Kit
- Mitochondrial & Genomic DNA Isolation Kits
- 5 Minutes DNA Ligation Kit
- 20 Minutes Gel Staining/Destaining Kit
- β-Galactosidase Staining Kit & Luciferase Reporter Assay Kit

Growth Factors and Cytokines

Monoclonal and Polyclonal Antibodies