



Antioxidant Assay Kit

Catalog Number KA1622

100 assays

Version: 04

Intended for research use only

www.abnova.com

Table of Contents

Introduction	3
Intended Use	3
Background	3
Principle of the Assay	3
General Information	4
Materials Supplied	4
Storage Instruction	4
Materials Required but Not Supplied	4
Precautions for Use	4
Assay Protocol	5
Sample Preparation	5
Assay Procedure	5
Data Analysis.....	6
Calculation of Results.....	6

Introduction

Intended Use

- Applications:
 - ✓ Direct Assays: serum, plasma, urine, saliva and other biological samples, food and beverages.
 - ✓ Drug Discovery/Pharmacology: effects of drugs on TAC.
- Key Features:
 - ✓ Sensitive and accurate. Use 20 μ L sample. Linear detection range from 1.5 to 1000 μ M Trolox equivalents.
 - ✓ Simple and high-throughput. The procedure involves addition of a single working reagent and incubation for 10 min. Can be readily automated as a high-throughput assay for thousands of samples per day.

Background

An ANTIOXIDANT is a molecule capable of slowing or preventing the oxidation of other molecules. Antioxidants protect the cells from damages by reactive oxygen species which are produced in oxidation reactions in the cell. Antioxidants can be small molecules such as glutathione, vitamins, or macromolecules such as catalase, glutathione peroxidase. As oxidative stress contributes to the development of many diseases including Alzheimer's disease, Parkinson's disease, diabetes, rheumatoid arthritis and neurodegeneration, the use of antioxidants in pharmacology is intensively studied. Antioxidants are also widely used as dietary supplements and in industry as preservatives in food, cosmetics, rubber and gasoline.

Simple, direct and high-throughput assays for total antioxidant capacity (TAC) find wide applications in research, food industry and drug discovery.

Principle of the Assay

The Antioxidant Assay Kit improved assay measures total antioxidant capacity in which Cu^{2+} is reduced by antioxidant to Cu^+ . The resulting Cu^+ specifically forms a colored complex with a dye reagent. The color intensity at 570 nm is proportional to TAC in the sample.

General Information

Materials Supplied

List of component

Component	Amount
Reagent A	12 mL
Reagent B	1 mL
Standard (50 mM Trolox)	100 µL

Storage Instruction

Store all components at -20 °C. Shelf life of six months after receipt.

Materials Required but Not Supplied

Pipetting devices, centrifuge tubes, clear flat-bottom uncoated 96-well plates, plate reader capable of reading optical density at 570 nm, homogenizer or sonicator etc.

Precautions for Use

Reagents are for research use only. Normal precautions for laboratory reagents should be exercised while using the reagents.

Assay Protocol

Sample Preparation

Samples should not contain any metal chelators (e.g. EDTA) and should be clear and free of any turbidity or particles. Liquid samples (e.g. non-hemolyzed serum, plasma) can be assayed directly. Cell lysate is prepared by homogenizing or sonicating cells in ice-cold 1 x PBS and centrifugation for 10 min at 14,000 rpm to pellet any debris. Use the clear supernatant for the assay. If not assayed immediately, freeze supernatant at -80°C (stable for 1 month).

Assay Procedure

1. Standards and Samples

Equilibrate all components to room temperature. Briefly centrifuge Reagent B and Standard before opening.

Mix 5 µL of the standard with 245 µL dH₂O (final 1 mM Trolox). Dilute standards as shown in the Table below. Transfer 20 µL standards into wells of a clear flat-bottom 96-well plate.

No	1 mM Trolox + H ₂ O	Vol (µL)	Trolox (µM)
1	100 µL + 0 µL	100	1000
2	60 µL + 40 µL	100	600
3	30 µL + 70 µL	100	300
4	0 µL + 100 µL	100	0

Transfer 20 µL of each sample into separate wells of the 96-well plate.

Note: for unknown samples, perform several dilutions to ensure that TAC is within the linear range of 1.5 to 1000 µM Trolox equivalents.

2. Assay

Prepare enough Working Reagent for Sample and Standard wells by mixing, for each assay well, 100 µL Reagent A and 8 µL Reagent B. Add 100 µL Working Reagent to all assay wells. Tap plate to mix. Incubate 10 min at room temperature.

3. Read OD_{570 nm} on a plate reader

Note: if calculated TAC is higher than 1000 µM Trolox equivalents, dilute sample in dH₂O and repeat assay. Multiply the results by the dilution factor.

Data Analysis

Calculation of Results

Subtract blank OD value (#4) from all standard and sample OD values. Plot the $\Delta OD_{570\text{ nm}}$ against standard concentrations and determine the slope of the standard curve. Calculate the Total Antioxidant Capacity (TAC) of Sample,

$$\text{TAC} = \frac{\text{OD}_{\text{SAMPLE}} - \text{OD}_{\text{BLANK}}}{\text{Slope } (\mu\text{M}^{-1})} \times n \text{ (}\mu\text{M Trolox Equivalent s)}$$

$\text{OD}_{\text{SAMPLE}}$ and OD_{BLANK} are the $\text{OD}_{570\text{ nm}}$ values of the sample and H_2O blank (standard #4). n is the sample dilution factor.

